

REF: P60108

FIBROBLAST MEDIUM

Product Type: Fibrobast Medium (w/o FBS)

Catalog Number: P60108 SF

国内流通品はFBSを含んでおりません。別途、FBSを10mL添加してご使用ください。

Product Description

Fibroblast Medium (FM) is a complete medium designed for optimal growth of normal human fibroblasts in vitro. It is a sterile, liquid medium which contains essential and non-essential amino acids, vitamins, organic and inorganic compounds, hormones, growth factors, trace minerals and a low concentration of fetal bovine serum (2%). The medium is HEPES and bicarbonate buffered and has a pH of 7.4 when equilibrated in an incubator with an atmosphere of 5% CO₂/95% air. The medium is formulated (quantitatively and qualitatively) to provide a defined and optimally balanced nutritional environment that selectively promotes proliferation and growth of normal human fibroblasts in vitro.

© Components

- 500 ml of Basal Medium
- 5 ml of Fibroblast Growth Supplement (FGS)
- 5 ml of penicillin/streptomycin solution (P/S solution)

別途、10mLのFBSを添加してください。

Prepare for use

Thaw FGS, FBS and P/S solution at 37°C. Gently tilt the FGS tube several times during thawing to help the contents dissolve. Make sure the contents of the supplement are completely dissolved into solution before adding to the medium. Rinse the bottle and tubes with 70% ethanol, and then wipe to remove excess. Remove the cap, being careful not to touch the interior threads with fingers. Add FGS. FBS and P/S solution into basal medium in a sterile field, mix well and then the reconstituted medium is ready for use. Since several components of this medium are lightlabile, it is recommended that the medium not be exposed to light for lengthy periods of time. If the medium is warmed prior to use, do not exceed 37°C. When stored in the dark at 4°C, the reconstituted medium is stable for one month.

© Caution

If handled improperly, some components of the medium may present a health hazard. Take appropriate precautions when handling it, including the wearing of protective clothing and eyewear. Dispose of properly.

-661-0978 兵庫県尼崎市久々知西町2丁目1-20 (お問合せ先) TEL:06-6435-9747 FAX:06-6435-9748 URL: http://www.kacnet.co.jp/ E-mail: shiyaku-info@kacnet.co.jp