



REF: P60116

MICROGLIAL CELL MEDIUM

Product Type:Microglial Cell MediumCatalog Number:P60116国内流通品はFBSを含んでおりません。非働化したFBSを25mL添加してご使用ください。

Microglia Medium (MM) is a complete medium designed for optimal growth of normal human brain-derived microglia in vitro. It is a sterile, liquid medium which contains essential and non-essential amino acids, vitamins, organic and inorganic compounds, hormones, growth factors, trace minerals and a 5% concentration of fetal bovine serum. The medium is bicarbonate buffered and has a pH of 7.4 when equilibrated in an incubator with an atmosphere of 5% CO₂/95% air. The medium is formulated (quantitatively and qualitatively) to provide a defined and optimally balanced nutritional environment that selectively promotes growth of normal human microglia in vitro.

\delta Components

- 500 ml of Basal Medium
- 5 ml of Microglia Growth Supplement (MCGS).
- 5 ml of penicillin/streptomycin solution (P/S solution)

別途、非働化FBS (HD-FBS) を25mL添加して ください。

🧔 Prepare for use

Thaw MGS, HD-FBS and P/S solution at 37°C. Gently tilt the MGS tube several times during thawing to help the contents dissolve. Make sure the contents of the supplement are completely dissolved into solution before adding to the medium. Rinse the bottle and tubes with 70% ethanol, and then wipe to remove excess. Remove the cap, being careful not to touch the interior threads with fingers. Add MGS, HD-FBS and P/S solution into basal medium in a sterile field, mix well and then the reconstituted medium is ready for use. Since several components of this medium are light-labile, it is recommended that the medium not be exposed to light for lengthy periods of time. If the medium is warmed prior to use, do not exceed 37°C. When stored in the dark at 4°C, the reconstituted medium is stable for one month.

S Product Use

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. Not approved for human or veterinary use, for application to humans or animals, or for use in vitro diagnostic or clinical procedures. 株式会社ケー・エー・シー

(ホスム) エン・シン 試薬事業部 〒661-0978 兵庫県尼崎市久々知西町2丁目1-20

(お問合せ先) TEL:06-6435-9747 FAX:06-6435-9748 URL:http://www.kacnet.co.jp/ E-mail:cs-info@kacnet.co.jpC

INNOVATIVE TECHNOLOGIES IN BIOLOGICAL SYSTEMS, S.L. Parque Tecnológico Bizkaia, Edf. 502, 1ª Planta | 48160 | Derio | Bizkaia Tel.: +34 944005355 | Fax: +34 946579925 innoprot@innoprot.com | www.innoprot.com